Compass

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60031632 Compass Data Management Software

PC system requirements:

Note that the below stated requirements are only for the Compass Application/Database and that requirements for the users operating system might be higher.

Compass Application

- Windows® 2000 Service Pack 4/Windows® Server 2003 Service Pack 1/Windows® XP Service Pack 2/Windows® Vista with SP 1
- Pentium III 600 MHz or better (1 GHz recommended)
- 192 MB RAM (512 MB recommended)
- 1025 MB available space on hard drive
- One Ethernet port for communication with the Analyzer

Compass Database

- Windows® 2000 Service Pack 4/Windows® Server 2003 Service Pack 1/Windows® XP Service Pack 2/Windows® Vista with SP 1
- .NET Framework 2.0¹⁾
- Microsoft SQL Server 2005 Express
- Pentium III 600 MHz or better (1 GHz recommended)
- 192 MB RAM (512 MB recommended)
- 1025 MB available space on hard drive
- One Ethernet port for communication with Compass Application

¹⁾Can be downloaded from Microsoft® website. .NET Framework 2.0 should be installed before Microsoft® SQL Server 2005 Express

Balance communication:

The PC should be connected to the balance via a serial cable (RS-232) according to the balance vendor's specification for PC communication (not included).

Listed commands for the following balance models are supported:

Sartorius A120S, A200S, A210P Models

GetWeight: ESC P CR LF
GetStableWeight: ESC P CR LF

Sartorius CP, GC and GP Models

GetWeight: ESC P CR LF
GetStableWeight: ESC P CR LF

Mettler Toledo Excellence Plus XP Models (XP105DR, XP205, XP205DR, XP204, XP504, XP504DR)

GetWeight: SI CR LF
GetStableWeight: S CR LF

LIMS communication

Export of result data to files in format *.txt, *.xml or *.csv for reading by LIMS

Import of sample data in file format *.xml created by LIMS

Required cables, not supplied in Compass package:

- Serial cable – (RS-232) connection to balance according to specification from balance vendor
- Network cable – Installation in network
- Crossover cable – Direct connection to Analyzer

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